

Wasp Plate Sealer

Outer cover removal procedure

1. Disconnect power and Pneumatic services to Wasp Unit (Ensure Heater block is COOL)
2. Remove foil reel from rear of Unit.
3. Remove the 4 outer cover bottom screws to free the main cover from the inner base unit
4. Remove 3 screws and remove left hand side inspection cover.



5. Push Heater block down (**ensure it is cool before touching it**)



6. Remove the screen Bezel Assembly



Wasp

Outer cover removal procedure

7. To remove display, Tilt display backwards in opening and very carefully withdraw from cover.



8. Disconnect all connections from screen
 - a. Screen Ribbon Cable.
 - b. Two Pin Fan Lead
 - c. Thermocouple (two pin black connector Pulls off) Green wire is + positive
 - d. Unplug the screen earth cable (green/yellow cable)



9. Remove display from cover.
10. Extreme care to be taken, do not separate the two Halves of touch screen!



Wasp

Outer cover removal procedure

11. Carefully lift cover from main chassis unit.



12. To replace cover, reverse this procedure.